

UV Imaging

Wiktor

July 19, 2010

Where In the room right of the coffee machine. Next to the OD machine/near the ice box.

Material needed

Quantity	What
	UV machine
	Gel, which you want to image.

Steps

1. Put the gel into the UVbox. Place it carefully to avoid bubbles close to the region we are interested in.
2. Close the door and turn on the UV light.
3. Computer: Open Kodak.M1.exe (Use guest account if nobody is logged in.)
4. "Capture GL 100" → "Preview" → "Capture".
5. We can adjust brightness or invert colors as desired.
6. "File" → "Export data" → "image" → save it on a memory stick.
7. Turn off the UV, and throw it away in the "white trash" under the gel bench.
8. We can pour the 0.5XTAE which is left in the box back into the bottle.

Warnings

- Do not touch computer keyboard with gloves!!
- Don't look at UV light! Don't turn UV on when box is not closed.