

LB agar medium recipe

Stefano Varricchio - EPFL iGEM 2010

July 15, 2010

Where On the wetbench at the right of the fume hood

		Quantity	What
Material needed	For 1 liter:	10 g	Bacto tryphone
		5 g	Yeast extract
		10 g	NaCl
		15 g	Agar
		1 L	diH ₂ O
		1	Magnetic stirrer

Steps

1. mix everything together and add the magnetic stirrer.
2. Sterilize by autoclaving.

Warnings

- Before doing a lot of bottles of medium try when they fit in the autoclave
 - Don't close the lid hermetically for autoclaving, the bottle will *explode*. Let air flow in and out the bottle.
 - Bacto yeast stinks a lot.
-