

iGEM TU/e 2014
Biomedical Engineering

Eindhoven University of Technology
Room: Ceres 0.04
Den Dolech 2, 5612 AZ Eindhoven
The Netherlands
Tel. no. +31 50 247 55 59
2014.igem.org/Team:TU_Eindhoven

Plasmid and gene digestion

Table of contents

Title Plasmid and gene digestion	1	Plasmid and gene digestion	3
--	----------	-----------------------------------	----------

1 Plasmid and gene digestion

Component	Quantity/mass/final concentration	Volume (μL)
H ₂ O	Fill up to 50 μL	..
10x buffer (Cut-Smart)	1 X	5
Plasmid DNA	5 μg (stock ... ng/ μL)	..
Restriction enzyme 1	20 U (20 U/ μL stock)	1
Restriction enzyme 2	20 U (20 U/ μL stock)	1
<i>Total</i>		50

- Mix well by pipetting up and down
- Incubate for 1 hour at 37°C, in PCR machine
- Heat inactivate for 20 min at 65°C, in PCR machine
- Store in fridge (-30°C)