

iGEM TU/e 2016
Biomedical Engineering

Eindhoven University of Technology
Room: Ceres 0.04
Den Dolech 2, 5612 AZ Eindhoven
The Netherlands
Tel. no. +31 50 247 55 59
2016.igem.org/Team:TU-Eindhoven

Transformation into NovaBlue

Table of contents

| | | | |
|------------------------------|----------|--|----------|
| Transformation into NovaBlue | 1 | Transformation into NovaBlue | 3 |
| | 1.1 | Materials | 3 |
| | 1.2 | Protocol | 3 |
| | 2 | Plating of the cells on agar plates | 3 |

1 Transformation into NovaBlue

Estimated bench time: 30 minutes

Estimated total time: 1.5 hour

Purpose: Amplification of the ligation product

It is essential to work sterile, thus disinfect your hands and work near a Bunsen Burner.

1.1 Materials

- Bucket with ice
- Bunsen Burner
- Eppendorf tubes with 20 mL competent NovaBlue cells
- Heat/shaking-block
- Incubator
- LB-agar plates supplemented with the correct antibiotic
- Ligation mixture
- Pipettes and tips
- SOC solution (Super optimal broth with catabolite repression)
- Water bath

1.2 Protocol

- Switch on the water bath and set temperature at 42 °C. Also turn on the heat/shaking-block and set up to 37 °C.
- Load a bucket with ice from the ice machine.
- Take the bacterial cells and SOC out of the -80 °C freezer. Transfer the cells directly to ice. Do not touch the bottom of the tube that contains the cells.
- Thaw the cells on ice for ~5 minutes.
- Add 1 µL of ligation mixture to 20 µL bacteria. Leave the bacteria on ice. Mix well. Make sure you work near the Bunsen burner flame.
- Incubate on ice for 5 min.
- Heat shock the cells for exactly 30 seconds at 42 °C.
- Return the cells directly to ice for 2 minutes.
- Add 80 µL of SOC solution (room temperature) to the bacteria. Do not return to ice!
- Incubate for 60 min at 37 °C and 300 rpm.
- Dry the LB-agar plates supplemented with the correct antibiotic in the incubator. Place the plates upside down (with the agar up) and slightly opened.

2 Plating of the cells on agar plates

Estimated bench time: 15 minutes

Estimated total time: 18 hour

Purpose: Amplification of the ligation product.

For more information, see our general Plating protocol.