

SOP Name: T4 Ligase

Author: Marcia Pryce

Source(s): Adapted from Exeter 2015

Ligase Materials:

COMPONENT 20 μ l REACTION

T4 DNA Ligase Buffer (10X)* 2 μ l

Vector DNA (4 kb) 50 ng (0.020 pmol)

Insert DNA (1 kb) 37.5 ng (0.060 pmol)

Nuclease-free water to 20 μ l

T4 DNA Ligase 1 μ l

Procedure:

1. Set up the following reaction in a microcentrifuge tube on ice. (T4 DNA Ligase should be added last. Note that the table shows a ligation using a molar ratio of 1:3 vector to insert for the indicated DNA sizes.) Use NEBioCalculator to calculate molar ratios. The T4 DNA Ligase Buffer should be thawed and resuspended at room temperature.
2. Gently mix the reaction by pipetting up and down and microfuge briefly.
3. For cohesive (sticky) ends, incubate at 16°C overnight or room temperature for 10 minutes.
4. For blunt ends or single base overhangs, incubate at 16°C overnight or room temperature for 2 hours (alternatively, high concentration T4 DNA Ligase can be used in a 10 minute ligation).
5. Heat inactivate at 65°C for 10 minutes.
6. Chill on ice and transform 1-5 μ l of the reaction into 50 μ l competent cells.