

Ni-NTA Spin Column Purification of 6xHis-Tagged Proteins under Native Conditions from *E. coli* Cell Lysates

The Ni-NTA Spin Kit (cat. no. 31314) and Ni-NTA Spin Columns (cat. no. 31014) can be stored at 2–8°C for up to 18 months.

For more information, please refer to the *Ni-NTA Spin Kit Handbook*, which can be found at www.qiagen.com/handbooks.

For technical assistance, please call toll-free 00800-22-44-6000, or find regional phone numbers at www.qiagen.com/contact.

Notes before starting

- Prepare bacterial cell pellet (stored at –20°C or –80°C).
 - Buffer compositions are provided in the appendix of the *Ni-NTA Spin Kit Handbook*.
 - Lysis requires Benzonase® nuclease (e.g., Novagen cat. no. 70664-3).
1. Resuspend a pellet derived from 5 ml cell culture volume in 630 µl Lysis Buffer (NPI-10). Add 70 µl Lysozyme Stock Solution (10 mg/ml) and add 3 Units/ml culture volume Benzonase Nuclease (i.e., for cell pellets from 5 ml cultures, add 15 Units Benzonase Nuclease).
 2. Incubate on ice for 15–30 min.
 3. Centrifuge lysate at 12,000 x g for 15–30 min at 4°C. Collect supernatant.
Note: Save 20 µl of the cleared lysate for SDS-PAGE analysis.
 4. Equilibrate the Ni-NTA spin column with 600 µl Buffer NPI-10. Centrifuge for 2 min at 890 x g (approx. 2900 rpm).
Note: The spin columns should be centrifuged with an open lid to ensure that the centrifugation step is completed under 2 min.

January 2011



5. Load up to 600 μ l of the cleared lysate containing the 6xHis-tagged protein onto the pre-equilibrated Ni-NTA spin column. Centrifuge for 5 min at 270 x g (approx. 1600 rpm), and collect the flow-through.

Note: To ensure efficient binding, do not exceed 270 x g when centrifuging Ni-NTA spin columns. The spin columns can be centrifuged with an open lid.

Note: Save the flow-through for analysis by SDS-PAGE.

6. Wash the Ni-NTA spin column twice with 600 μ l Buffer NPI-20. Centrifuge for 2 min at 890 x g (approx. 2900 rpm).

Note: Save the flow-through for analysis by SDS-PAGE.

7. Elute the protein twice with 300 μ l Buffer NPI-500. Centrifuge for 2 min at 890 x g (approx. 2900 rpm), and collect the eluate.

Note: Most of the 6xHis-tagged protein (>80%) should elute in the first 300 μ l eluate. The remainder will elute in the second 300 μ l.

For up-to-date licensing information and product-specific disclaimers, see the respective QIAGEN kit handbook or user manual.

Trademarks: QIAGEN® (QIAGEN Group); Benzonase® (Merck KGaA).
1067581 01/2011 © 2011 QIAGEN, all rights reserved.

