



## SOC-medium

- ◆ Add the following components for 900 mL of distilled H<sub>2</sub>O
  - ◇ 20 g Trypton
  - ◇ 5 g Bacto Yeast Extract
  - ◇ 2 mL of 5 M NaCl
  - ◇ 2.5 mL of 1 M KCL
  - ◇ 10 mL of 1 M MgCl<sub>2</sub>
  - ◇ 10 mL of 1 M MgSO<sub>4</sub>
  - ◇ 20 mL of 1 M glucose

From: [iGEM Bielefeld-CeBiTec 2014](#)