

07. (July) 2019

Project: iGEM_Munich2019 Shared Project

Authors: Theresa Keil

WEDNESDAY, 17/7/2019

- SDS-Page: samples were prepared on 16.07.19
- run: 200 V 35 min
- Western Blot protocol: put chamber on ice to prevent overheating, run 100 V, 1 h
- washed membrane with water after run, continued with ponceaus staining
- Blocking in 5 % milk in TBS for 1 h at RT
- Incubation in 1st AB mouse-alpha-CD63 (2µg/ml) in 5 % milk in TBS over night at 4 °C