

Jul 10, 2019

## 14 SDS-PAGE

TJUSLS China<sup>1</sup><sup>1</sup>Tianjin University

1

Works for me

dx.doi.org/10.17504/protocols.io.49qgz5w

TJUSLS China  
Tianjin University

## GUIDELINES

SDS-PAGE Separation of gel samples	Scope of application
6% sample gel	50-150kD
8% sample gel	30-90kD
10% sample gel	20-80kD
12% sample gel	12-60kD
15% sample gel	10-40kD

## MATERIALS

NAME	CATALOG #	VENDOR
TEMED	1610801	Bio-rad Laboratories
40g Acrylamide/ Bisacrylamide (29:1); Premixed Powder	786-506	G-Biosciences
Coomassie brilliant blue G-250	CB0038.SIZE.100g	Bio Basic Inc.
SDS, 10%(w/v) solution	SD8118.SIZE.100ml	Bio Basic Inc.
Tris pH 6.8		
10% Ammonium persulfate	/	
1M Tris(PH=8.8)	/	

## SAFETY WARNINGS

Be careful of 30% Acr-Bis M and TEMED owing to their potential harm.

## BEFORE STARTING

In general, 15% sample gel is used, so here is only the composition of 15% sample gel.


**Composition of the 15% separated of the gel sample(4mL):**

Composition	volume/mL
Distilled water	0.4
30% Acr-Bis 29:1	2
1M Tris pH=8.8	1.5
10% SDS	0.04
10% Ammonium persulfate	0.04
TEMED	0.002

**Composition of concentrated sample gel(1mL):**

Composition	volume/mL

Distilled water	0.7
30% Acr-Bis	0.165
1M Tris pH=6.8	0.125
10% SDS	0.01
10% Ammonium persulfate	0.01
TEMED	0.001

- 1 Mount the gels in the vertical electrophoresis apparatus.
- 2 Load samples under the cathode buffer. Apply 10 $\mu$ L sample volumes to 0.7  $\times$  5 mm sample wells.  
 10  $\mu$ l
- 3 Set running conditions appropriate to your type of gel. We usually set 200V to run at the beginning, after 10 minutes change the voltage into 300V for another 30 minutes.
- 4 Protein can be visualized directly in the gel by Coomassie staining or silver staining. We usually choose Coomassie staining, then transfer the gel to the decolorizing solution and remove the dye solution.
- 5 Recording with gel imager.



This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited