05. (May) 2019

Project: iGEM_Munich2019 Shared Project

Authors: Sarah Brajkovic

MONDAY, 6/5/2019

Cell splitting:

HEK 293 T cells were splitted according to cell splitting protocol.

Measurements from cell counting: 3.01 x10^6 cells/mL

Mixed 170 µl of cell suspension to 12 ml Advanced DMEM with FBS and put 1 ml in each well

Seeding:

HEK 293 T cells were seeded onto a 12-well plate.

12 wells were seeded with each well containing 0.1 x 10^6 cells/mL

Incubation at 37°C, 5% CO2

file:///tmp/tmpRK1GNI.html