

08. (August) 2019

Project: iGEM_Munich2019 Shared Project

Authors: Theresa Keil

WEDNESDAY, 14/8/2019

cells from VLP team

washed 3 times with PBS, spin down and discard supernatant

lysis with 1 % triton 60 °C, 10 min

DNase treatment

continued with trizol treatment (incubated with trizol 10 min)

tried to increase 260/230 ratio by washing twice with ethanol

sample 1: 14.7 ng/μl , 260/280=1.56, 260/230=0.51

sample 2: 44.9 ng/μl, 260/280=1.61, 260/230=0.27