## 05. (May) 2019

Project: iGEM\_Munich2019 Shared Project

Authors: Sarah Brajkovic

MONDAY, 13/5/2019

## Cell splitting:

HEK 293 T cells were splitted according to cell splitting protocol.

Measurements from cell counting: 3 x10^6 cells/mL

Mixed 400 µl of cell suspension to 12 ml Advanced DMEM with FBS and put 1 ml in each well

## Seeding:

HEK 293 T cells were seeded onto a 12-well plate.

12 wells were seeded with each well containing 0.1 x 10^6 cells/mL

Incubation at 37°C, 5% CO2

file://tmp/tmpgrAtAV.html