



Jul 08, 2019

03 Ligation

TJUSLS China¹¹Tianjin University

1

Works for me

dx.doi.org/10.17504/protocols.io.49iqz4e

TJUSLS China
Tianjin University

MATERIALS

NAME ▾

CATALOG # ▾

VENDOR ▾

ddH₂O

Water bath or heat block at 37°C

5×T4 DNA Ligase Buffer

/

T4 DNA Ligase

/

- Mix together:

5×T4 DNA Ligase Buffer	2μl
T4 DNA Ligase	0.5–1μl
Vector	as required
Insert	as required
ddH ₂ O	Add to 10μl

- Heat in water bath or heat block at 22°C for 30min.

22 °C

00:30:00



This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited