



Oct 06, 2019

17 Inhibition Kinetics Measurement V.2

TJUSLS China¹¹Tianjin University

1

Works for me

dx.doi.org/10.17504/protocols.io.7y9hpb6

TJUSLS China
Tianjin University

MATERIALS

NAME ▾	CATALOG # ▾	VENDOR ▾
DMSO	GB-D-360	P212121
Scale		
General 96-well plates (Black)	/	
Infinite M1000 Pro Automatic Microplate Reader	/	
Multi-channel adjustable pipette	/	
Fluorescent Probe(CDC-1)	/	
Target Enzyme(beta-lactamase)	/	
Ultrasonic Cleaner		

BEFORE STARTING

Reference

[1] Cer RZ, Mudunuri U, Stephens R, Lebeda FJ. IC50-to-Ki: a web-based tool for converting IC50 to Ki values for inhibitors of enzyme activity and ligand binding. Nucleic Acids Res. 2009;37(Web Server issue):W441–W445. doi:10.1093/nar/gkp253

IC50 in vitro test

- 1 Soak the 96-well plates in 75% ethanol and put the container in ultrasonic cleaner for 30min to 1 hour, then use ddH₂O to wash these plates several times. Put these clean plates in drying oven at 55°C.
- 2 Dilute inhibitors using 100%DMSO to 8mM (a.k.a. 80μM in reaction system) and dilute in 2-time gradients for 12 concentrations. Dilute the enzyme using its buffer to standard concentration (same as the determination of Km).
- 3 Add the reaction system into 96-well plates.
Pipet 94μL protein solution and 1μL inhibitor solution(diluted as 2-time gradients, at least 8 different concentrations), incubated at room temperature for 5 mins, then add substrate CDC-1 5μL.
- 4 Set controls.
Negative control: 94μL protein solution+1μL DMSO, 5μL substrate
Blank control: 94μL protein buffer+1μL DMSO, 5μL substrate
And there are 3 parallel holes for each system.

- 5 Set up the program in Infinite M1000 Pro Automatic Microplate Reader.
Shake for 10 sec at 654 rpm
Kinetic Cycle (to read fluorescent intensity each cycle)
Fluorescent measure, 20 cycle, 30sec for each cycle
- 6 Put the plate in Microplate reader, and click Start button.
- 7 When the facility ends testing, save data and import the data of 0-200s into GraphPad Prism Software. Use "nonlinear fit" – "straight line" and regulate the number of data to fit R^2 .
- 8 Move baseline of blank control. Calculate $I_r = (1 - V_r/V_0) \times 100\%$. Take average I_r of each inhibitors' concentrations as Y value, and take $\log[I]$ as X value. Then use "nonlinear fit" – "log(inhibitor) vs. normalized response – Variable slope" to fit IC_{50} curve, and its value would be calculated automatically.

Judging of inhibition mechanism

9 reversible/irreversible inhibition

[Set up gradient concentrations of protein\(revolves around the value in screening system\)](#)

As well as gradient concentrations of inhibitor(revolves around the value of IC_{50})

[Draw the plot of \$V_0-\[E\]\$ to see the movement of straight line.](#)

If they are parallel with each other, the mechanism is irreversible.

If they are all through the origin, the mechanism is reversible.

10 competitive/noncompetitive inhibition

Set up gradient concentrations of substrate(revolves around the value in screening system)

As well as gradient concentrations of inhibitor(revolves around the value of IC_{50}).

Draw the plot of $1/V_0 - 1/[S]$ to see the movement of straight line (Lineweaver-Burk plot).

If they gather at Y axis, the mechanism is competitive.

If they pass the same point in X axis, the mechanism is noncompetitive.

If they gather at another point in this dimension, it would be mixed type competition.(uncompetitive inhibition)

11 K_i

After decision of this inhibitor's mechanism, a constant K_i can be calculated via different equations.^[1]

For competitive inhibition,

$$K_i = \frac{IC_{50}}{(S/K_m + 1)} \begin{cases} \text{if } S = K_m, & K_i = IC_{50}/2 \\ \text{if } S \gg K_m, & K_i \ll IC_{50} \\ \text{if } S \ll K_m & K_i \cong IC_{50} \end{cases}$$

For noncompetitive inhibition,

$$K_i = IC_{50} \text{ when } S = K_m \text{ or } S \gg K_m \text{ or } S \ll K_m$$

For uncompetitive inhibition,

$$K_i = \frac{IC_{50}}{(K_m/S + 1)} \begin{cases} \text{if } S = K_m, & K_i = IC_{50}/2 \\ \text{if } S \gg K_m, & K_i \cong IC_{50} \\ \text{if } S \ll K_m & K_i \ll IC_{50} \end{cases}$$

Also, we use a website tool (https://bioinfo-abcc.ncicrf.gov/IC50_Ki_Converter/index.php) to calculate its value automatically.



This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited