

08. (August) 2019

Project: iGEM_Munich2019 Shared Project

Authors: Johanna Wallner

TUESDAY, 13/8/2019

Johanna

cell culture: transfection for Purification and qPCR

- 550 µL medium were exchanged before transfection
- 2 24-well plates seeded the 12th of August

Transfection scheme - qPCR and Purification 13/08/19 (ng per well)


	Condition	V8	V10	V11	V14	V15	V27	V28	V30
1	2	200 ng	-	-	-	-	-	-	300 ng
2	4	200 ng	-	100 ng	100 ng	-	100 ng	-	-
3	6	200 ng	100 ng	-	100 ng	-	100 ng	-	-
4	7	200 ng	100 ng	-	-	100 ng	100 ng	-	-
5	9	200 ng	-	100 ng	-	100 ng	100 ng	-	-
6	4H	200 ng	-	100 ng	100 ng	-	-	100 ng	-
7	V14	-	-	-	500 ng	-	-	-	-

- Transfection:

Transfectionmix 24-well plate 13/0...

	A	B
1	DNA per well	500 ng
2	P3000 Reagent per well	1 µL
3	Lipofectamine 3000 reagent per well	0.75 µL
4	OptiMEM per well	2 x 25 µL

- Transfect cells according to the following table. Use the indicated volume of DNA and P3000™ Reagent with each of the two volumes of Lipofectamine™ 3000 (when performing optimization). Each reaction mix volume is for one well and accounts for pipetting variations. Scale volumes proportionally for additional wells.

 image.png

Timeline	Steps	Procedure Details (Two Reaction Optimization)			
		Condition 1	Condition 2	Condition 3	Condition 4
Day 0	1. Seed cells to be 70-80% confluent at transfection	Adherent cells	1.4 x 10 ⁶	0.5-2 x 10 ⁶	0.25-1 x 10 ⁶
	2. Dilute Lipofectamine [®] 3000 Reagent in Opti-MEM [®] Medium (2 tubes) - Mix well	Opti-MEM [®] Medium	5 µL x 2	25 µL x 2	125 µL x 2
		Lipofectamine [®] 3000 Reagent	0.15 and 0.3 µL	0.75 and 1.5 µL	3.75 and 7.5 µL
Day 1	3. Prepare master mix of DNA by diluting DNA in Opti-MEM [®] Medium, then add P3000 [®] Reagent - Mix well	Opti-MEM [®] Medium	10 µL	50 µL	250 µL
		DNA (0.5-5 µg/µL)	0.2 µg	1 µg	5 µg
		P3000 [®] Reagent (2 µL/µg DNA)	0.4 µL	2 µL	10 µL
Day 1	4. Add Diluted DNA to each tube of Diluted Lipofectamine [®] 3000 Reagent (1:1 ratio)	Diluted DNA (with P3000 [®] Reagent)	5 µL	25 µL	125 µL
		Diluted Lipofectamine [®] 3000 Reagent	5 µL	25 µL	125 µL
	5. Incubate	Incubate for 10-15 minutes at room temperature.			
Day 2-4	6. Add DNA-lipid complex to cells	Condition 1 (per well)	Condition 2 (per well)	Condition 3 (per well)	Condition 4 (per well)
		DNA-lipid complex	10 µL	50 µL	250 µL
		DNA amount	100 ng	500 ng	2500 ng
Day 2-4	7. Visualize/analyze transfected cells	P3000 [®] Reagent	0.2 µL	1 µL	5 µL
		Lipofectamine [®] 3000 Reagent used	0.15 and 0.3 µL	0.75 and 1.5 µL	3.75 and 7.5 µL
		Incubate cells for 2-4 days at 37°C. Then, analyze transfected cells.			

- transfection finished at 10:00

Alejandro

cell culture: transfection for HiBit Assay in 96-well plates


- cells were at 50-60 % confluence (rather 50 % after medium exchange)
- 100 µL medium over the cells were exchanged before transfectin
- transfection was finished at 11:30
- 8 replicates

Transfection scheme - HiBit 96-well plate 13/08/19 (ng per well)								
	Condition	V8	V10	V11	V14	V15	V17	V30
1	1 -	-	-	-	-	-	-	100 ng
2	2 40 ng	-	-	-	-	-	20 ng	40 ng
3	4 40 ng	-	-	20 ng	20 ng	-	20 ng	-
4	6 40 ng	-	20 ng	-	20 ng	-	20 ng	-
5	7 40 ng	-	20 ng	-	-	20 ng	20 ng	-
6	9 40 ng	-	-	20 ng	-	20 ng	20 ng	-

- Transfection:

Transfectionmix 96-well plate 13/0...			A	B
1	DNA per well	100 ng		
2	P3000 Reagent per well	0.2 µL		
3	Lipofectamine 3000 reagent per well	0.15 µL		
4	OptiMEM per well	2 x 5 µL		

- Transfect cells according to the following table. Use the indicated volume of DNA and P3000[™] Reagent with each of the two volumes of Lipofectamine[™] 3000 (when performing optimization). Each reaction mix volume is for one well and accounts for pipetting variations. Scale volumes proportionally for additional wells.

 image.png

Timeline		Steps	Procedure Details (Two Reaction Optimization)																						
Day 1	1	Seed cells to be 70-90% confluent at transfection	<table><thead><tr><th>Component</th><th>24-well</th><th>96-well</th><th>48-well</th></tr></thead><tbody><tr><td>Adherent cells</td><td>1-4 x 10⁵</td><td>0.5-2 x 10⁵</td><td>0.25-1 x 10⁵</td></tr></tbody></table>				Component	24-well	96-well	48-well	Adherent cells	1-4 x 10 ⁵	0.5-2 x 10 ⁵	0.25-1 x 10 ⁵											
Component	24-well	96-well	48-well																						
Adherent cells	1-4 x 10 ⁵	0.5-2 x 10 ⁵	0.25-1 x 10 ⁵																						
2	Dilute Lipofectamine™ 2000 Reagent in Opti-MEM Medium (2 tubes) - Mix well	<table><tbody><tr><td>Opti-MEM® Medium</td><td>5 µL x 2</td><td>25 µL x 2</td><td>125 µL x 2</td></tr><tr><td>Lipofectamine™ 2000 Reagent</td><td>0.15 and 0.3 µL</td><td>0.75 and 1.5 µL</td><td>3.75 and 7.5 µL</td></tr></tbody></table>				Opti-MEM® Medium	5 µL x 2	25 µL x 2	125 µL x 2	Lipofectamine™ 2000 Reagent	0.15 and 0.3 µL	0.75 and 1.5 µL	3.75 and 7.5 µL												
Opti-MEM® Medium	5 µL x 2	25 µL x 2	125 µL x 2																						
Lipofectamine™ 2000 Reagent	0.15 and 0.3 µL	0.75 and 1.5 µL	3.75 and 7.5 µL																						
3	Prepare master mix of DNA by diluting DNA in Opti-MEM® Medium, then add P3000® Reagent - Mix well	<table><tbody><tr><td>Opti-MEM® Medium</td><td>10 µL</td><td>50 µL</td><td>250 µL</td></tr><tr><td>DNA (0.5-5 µg/µL)</td><td>0.2 µg</td><td>1 µg</td><td>5 µg</td></tr><tr><td>P3000® Reagent (2 µL/µg DNA)</td><td>0.4 µL</td><td>2 µL</td><td>10 µL</td></tr></tbody></table>				Opti-MEM® Medium	10 µL	50 µL	250 µL	DNA (0.5-5 µg/µL)	0.2 µg	1 µg	5 µg	P3000® Reagent (2 µL/µg DNA)	0.4 µL	2 µL	10 µL								
Opti-MEM® Medium	10 µL	50 µL	250 µL																						
DNA (0.5-5 µg/µL)	0.2 µg	1 µg	5 µg																						
P3000® Reagent (2 µL/µg DNA)	0.4 µL	2 µL	10 µL																						
4	Add diluted DNA to each tube of Diluted Lipofectamine™ 2000 Reagent (1:1 ratio)	<table><tbody><tr><td>Diluted DNA (with P3000® Reagent)</td><td>5 µL</td><td>25 µL</td><td>125 µL</td></tr><tr><td>Diluted Lipofectamine™ 2000 Reagent</td><td>5 µL</td><td>25 µL</td><td>125 µL</td></tr></tbody></table>				Diluted DNA (with P3000® Reagent)	5 µL	25 µL	125 µL	Diluted Lipofectamine™ 2000 Reagent	5 µL	25 µL	125 µL												
Diluted DNA (with P3000® Reagent)	5 µL	25 µL	125 µL																						
Diluted Lipofectamine™ 2000 Reagent	5 µL	25 µL	125 µL																						
5	Incubate	Incubate for 10-15 minutes at room temperature.																							
6	Add DNA-lipid complex to cells	<table><thead><tr><th>Component (per well)</th><th>96-well</th><th>24-well</th><th>6-well</th></tr></thead><tbody><tr><td>DNA-lipid complex</td><td>10 µL</td><td>50 µL</td><td>250 µL</td></tr><tr><td>DNA amount</td><td>100 ng</td><td>500 ng</td><td>2500 ng</td></tr><tr><td>P3000® Reagent</td><td>0.2 µL</td><td>1 µL</td><td>5 µL</td></tr><tr><td>Lipofectamine™ 2000 Reagent used</td><td>0.15 and 0.3 µL</td><td>0.75 and 1.5 µL</td><td>3.75 and 7.5 µL</td></tr></tbody></table>				Component (per well)	96-well	24-well	6-well	DNA-lipid complex	10 µL	50 µL	250 µL	DNA amount	100 ng	500 ng	2500 ng	P3000® Reagent	0.2 µL	1 µL	5 µL	Lipofectamine™ 2000 Reagent used	0.15 and 0.3 µL	0.75 and 1.5 µL	3.75 and 7.5 µL
Component (per well)	96-well	24-well	6-well																						
DNA-lipid complex	10 µL	50 µL	250 µL																						
DNA amount	100 ng	500 ng	2500 ng																						
P3000® Reagent	0.2 µL	1 µL	5 µL																						
Lipofectamine™ 2000 Reagent used	0.15 and 0.3 µL	0.75 and 1.5 µL	3.75 and 7.5 µL																						
Day 2	7	Visualize/measure transfected cells	Incubate cells for 2-4 days at 37°C. Then, analyze transfected cells.																						