

# 09. (September) 2019

---

**Project:** iGEM\_Munich2019 Shared Project

**Authors:** Theresa Keil

FRIDAY, 20/9/2019

---

- Receiving Min6 exosomes (starting volume: 1mL) + Anja 72h (starting volume: 290µl)
- Put Anajs 48h and 72h and Min6 in Exosome isolation kit and incubate in 4°C overnight
- RNA isolation out of HEK exosomes from 19.09. with Exosomal RNA and protein isolation kit; the manufacturer's protocol was slightly modified: the empty centrifugation after the washing steps was extended to 4 min, 13.000 rpm, RT; finally, the samples were eluted in 40µl 95°C hot water