NucleoSpin® Plasmid EasyPure

I) Preparations

- Add total content of Liquid RNase A to Buffer A1, mix thoroughly, and store at 2–8 °C.
- Add indicated volume of 96–100 % ethanol to Buffer AQ (Concentrate).
- Before each preparation, check Buffer A2 for precipitated SDS. Dissolve any precipitate for several minutes at 30–40 °C and store buffer at room temperature.

II) Culture volume

Cells can be grown in LB or rich media (e.g., TB, CircleGrow) but make sure not to exceed the recommended culture volumes (see table below). Using too much cell mass will exhaust the lysis capacity of the buffer system and result in poor or even no yield. If nevertheless more cells have to be processed (e.g., for low copy-number plasmids), increase lysis buffers A1, A2, and A3 proportionally.

OD _{600 nm}	1	2	3	4	5	6
Max. volume	15 mL	8 mL	5 mL	4 mL	3 mL	2 mL

For further information, please download the **NucleoSpin® Plasmid EasyPure** user manual at the relating webpage www.mn-net.com/plasmid

III) Protocol overview

1. Pellet cells into a microcentrifuge tube and discard supernatant.
2. Add **150 µL Buffer A1** and vortex to resuspend cells completely.
3. Add **250 µL Buffer A2** and invert 5 times – **DO NOT VORTEX!**
4. Incubate up to **2 min** at **room temperature** to lyse cells.
5. Add **350 µL Buffer A3** and invert **until lysate has turned colorless**.
6. Centrifuge for **3 min** at **full speed** (> 12,000 x g) to pellet precipitate.
7. Put a NucleoSpin® Plasmid EasyPure Column into a Collection Tube (2 mL).
8. Load clear supernatant onto the spin column.
9. Centrifuge for **30 s** at **1,000–2,000 x g** and discard flow-through.
10. Add **450 µL Buffer AQ** to the spin column.
11. Centrifuge for **1 min** at **full speed** (> 12,000 x g) and discard collection tube.
REPEAT IF SPIN COLUMN IS STILL WET!
12. Put spin column into a new 1.5 mL microcentrifuge tube (not provided).
13. Add **50 µL Buffer AE** onto the membrane and **incubate for 1 min***.
14. Centrifuge for **1 min** at **full speed** (> 12,000 x g).

* For larger constructs or large amounts of plasmid, elution efficiency can be increased by incubating spin cup with hot buffer at 70 °C, increasing the elution volume, or eluting several times with fresh buffer.

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V) Product use restriction/warranty

NucleoSpin® Plasmid EasyPure is developed, designed, and sold for research purpose only. We offer a one year warranty from date of delivery that our products will conform to applicable specifications set forth in the product specifications if not sold to persons set forth in § 13 of German Civil Code. In such a case, the provisions of the German Civil Code shall be valid. This warranty does not cover defects in or damage to the products which are due to improper installation or maintenance, misuse, neglect or any use other than ordinary commercial application.

The seller shall not be liable for the goods being fit for a particular purpose unless otherwise agreed upon, to which the buyer intends to put them. This warranty is strictly exclusive. MACHERY-NAGEL MAKES no other warranty of any kind whatsoever, and SPECIFICALLY DISCLAIMS AND EXCLUDES ALL OTHER WARRANTIES OF ANY KIND OR NATURE WHATSOEVER; DIRECTLY OR INDIRECTLY; EXPRESS OR IMPLIED; INCLUDING; WITHOUT LIMITATION; AS TO THE SUITABILITY; REPRODUCIBILITY; DURABILITY; FITNESS FOR A PARTICULAR PURPOSE OR USE; MERCHANTABILITY; CONDITION; OR ANY OTHER MATTER WITH RESPECT TO MACHERY-NAGEL PRODUCTS OR THE SALE OF MACHERY-NAGEL PRODUCTS. If not superseded by these terms the current terms and conditions for the sale of goods (export version) are additionally valid.

VI) Ordering information

Product	Pack of	REF
NucleoSpin® Plasmid EasyPure	10 / 50 / 250 preps	740727.10 / .50 / .250
NucleoSpin® Plasmid Buffer Set (for the isolation of low copy plasmids)	1	740953
Liquid RNase A	250 mg	740397

MACHERY-NAGEL

www.mn-net.com